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Qingdao ChuangMeng Instrument Co., Ltd.

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Frequency Conversion HighSpeed Mixers

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©Copyright owned by Qingdao ChuangMeng Instrument Co., Ltd

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Carefully read this User Manual to learn how to install and use the product correctly. After reading, properly keep the User Manual as a reference for future maintenance and repair.

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Address: No. 7 Wenzhou Road, City Northern District, Qingdao City, China

Production base: No. 3 Xinghai Road, Liuting Street, Chengyang District, Qingdao

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The variable frequency high-speed mixer produced by Chuangmeng Instrument adopts advanced variable frequency technology and microcomputer technology. It has functions such as timing, fixed speed, single axis and multi axis simultaneous operation. The mixing blades are made according to API standards. It has the characteristics of simple operation, high starting torque, stable speed, reliability, and low noise. It is specialized experimental equipment for petroleum, exploration, chemical and other industries. Mainly used for high-speed mixing of various drilling fluids, chemical reagents, etc. for petroleum drilling, it is an ideal high-speed mixing equipment.

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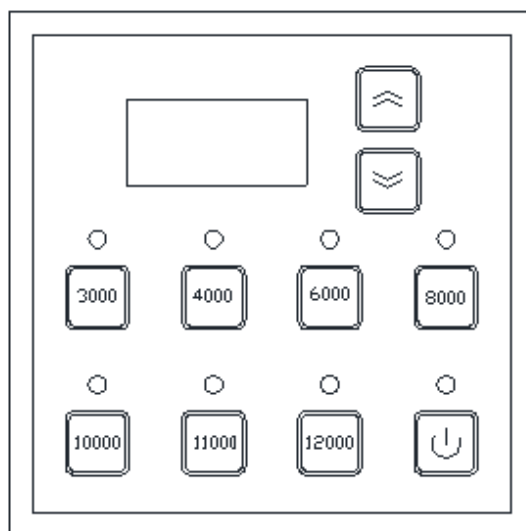
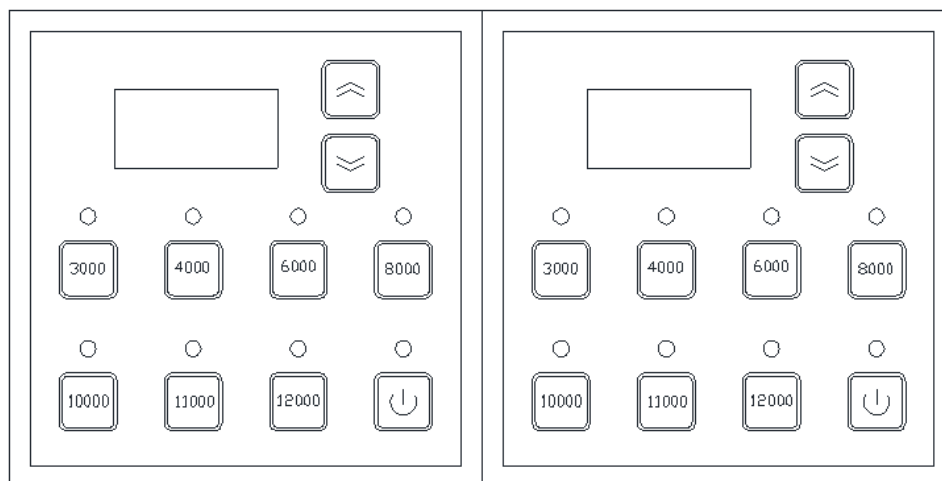
Schematic diagram of control box for 1613 high-speed mixer

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Schematic diagram of control box for 1614 high-speed mixer

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1. Connect the instrument to the corresponding voltage and power supply.

Note: the four axis mixer has two output lines, which should be connected to their respective mainframe respectively.

2. Pour the sample into the liquid cup. In order to prevent overflow when sample is stirred, the injection amount of sample is generally not more than 2/3.

3. Place the top of the liquid cup on the opening of the cup assembly, and place the cup on the lower support seat. The liquid cup is held tightly by the pressing cup assembly and the lower supporting seat.

4. When selecting the operation, select the gear according to the tachometer and set the mixing time. Finally, press the gear key to start the operation of the instrument, and the gear number flashes during the operation.
5. Select stop, (a) press the gear knob. (b) set the time until the instrument stops automatically. Note: gear position and timing can only be adjusted when the instrument is stopped.
6. After stirring, hold the cup tightly and lift it forcefully. Remove the cup from the lower support.

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- 1、清洗各部件并干燥待用，仪器置于干燥环境中。
 - 2、移动或保养仪器时。要轻拿、轻放，以免造成部件变形影响精度和使用。
 - 3、搅拌轴不得碰撞，以免弯曲变形，损坏仪器。
 - 4、电动机和调速器应经常保持清洁干燥不能受潮，要定期检查，严禁将腐蚀性化学药品与本机同存放。
 - 5、搅拌叶片若重量低于 5.1 克需更换。其原始重量约为 5.5 克。
1. Clean the components and dry them up. The instruments are placed in a dry environment.
 2. When moving, repairing or maintaining instruments. It is necessary to take lightly and put it lightly so as not to cause deformation of parts and affect accuracy and use.
 3. The stirring shaft must not collide, so as to avoid bending deformation

and damage the instrument.

4. Motor and speed regulator should be kept clean, dry and not damp. Regular inspection should be carried out. It is strictly forbidden to store corrosive chemicals with the machine.

5. If the weight of the stirring blade is less than 5.1g, it needs to be replaced. Its original weight is about 5.5g.

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故障 Fault	原因 Reason	维修方法 Maintenance method
电机噪音增大 Motor noise increase	电机损坏 Motor damage	更换电机
电机不转动 Motor does not rotate	①电机损坏 1. Motor damage ②线路板出现故障 2. Circuit board failure.	①更换电机 ②更换线路板 1. Replace the motor 2. Replacing circuit boards
只有部分转速 Only partial rotation speed	“转速转换按钮”损坏 "Speed conversion button" damage	更换转速转换按钮 Change the speed conversion button.
搅拌叶片不转 Stirring blades do not turn	紧固螺钉松动 Loosening of fastening screws	将紧固螺钉拧紧 Tighten fastening screws

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Production address: No. 3 Xinghai Road, Liuting Street, Chengyang District, Qingdao

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结论: Conclusion:

经检验, 青岛创梦仪器有限公司生产的产品符合上述标准的要求。准予出厂。
After inspection, Qingdao Chuangmeng Instrument Co., Ltd The products produced meet the requirements of the above standards. Approved for delivery.

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Qingdao Chuangmeng Instrument Co., Ltd. Packing list
Manufacturing enterprise: Qingdao Chuangmeng Instrument Co.,Ltd.